

Antigen-Specific T Cells for Assessing Immune Modulating Therapeutics

RoukenBio

BIOIVT
ELEVATING SCIENCE®

Matt Freer, Hannah Findlay, Agapitos Patakas, Robert Benson, Emilio Cosimo, Adele Hannigan

RoukenBio, Rouken Discovery Centre, Braidhurst Industrial Estate, Motherwell, ML1 3ST, UK



Introduction


Antigen-specific T cells play a key role in the adaptive immune system, as they can identify and eliminate cells presenting their specific antigen. Their targeted nature makes them valuable for testing immune-modulating therapies designed to either boost or suppress immune activity in diseases like cancer and autoimmunity. A major obstacle in evaluating antigen-specific T cell responses is having readily available cells to study the impact of immune modulators in real time. To address this, we have developed well-characterised transgenic and endogenous antigen-specific 'Thaw & Go' T cell banks that facilitate the evaluation of new immune modulators.

Generation of TCR-T Cells

Donors were selected based on HLA genotype compatibility with the antigen of interest, ensuring TCR-T cells will recognise their specific antigen when presented by the specific HLA genotype.


CD4+ or CD8+ T cells were isolated from cryopreserved PBMC (supplied by BioIVT, LLC) and transduced with a lentiviral vector encoding the TCR construct. An upstream tag enabled selection of successfully transduced cells. The cells were expanded using a proprietary protocol, and then characterised and cryopreserved.

Characterisation of the cells included transgenic TCR expression, and T cell exhaustion and activation markers.

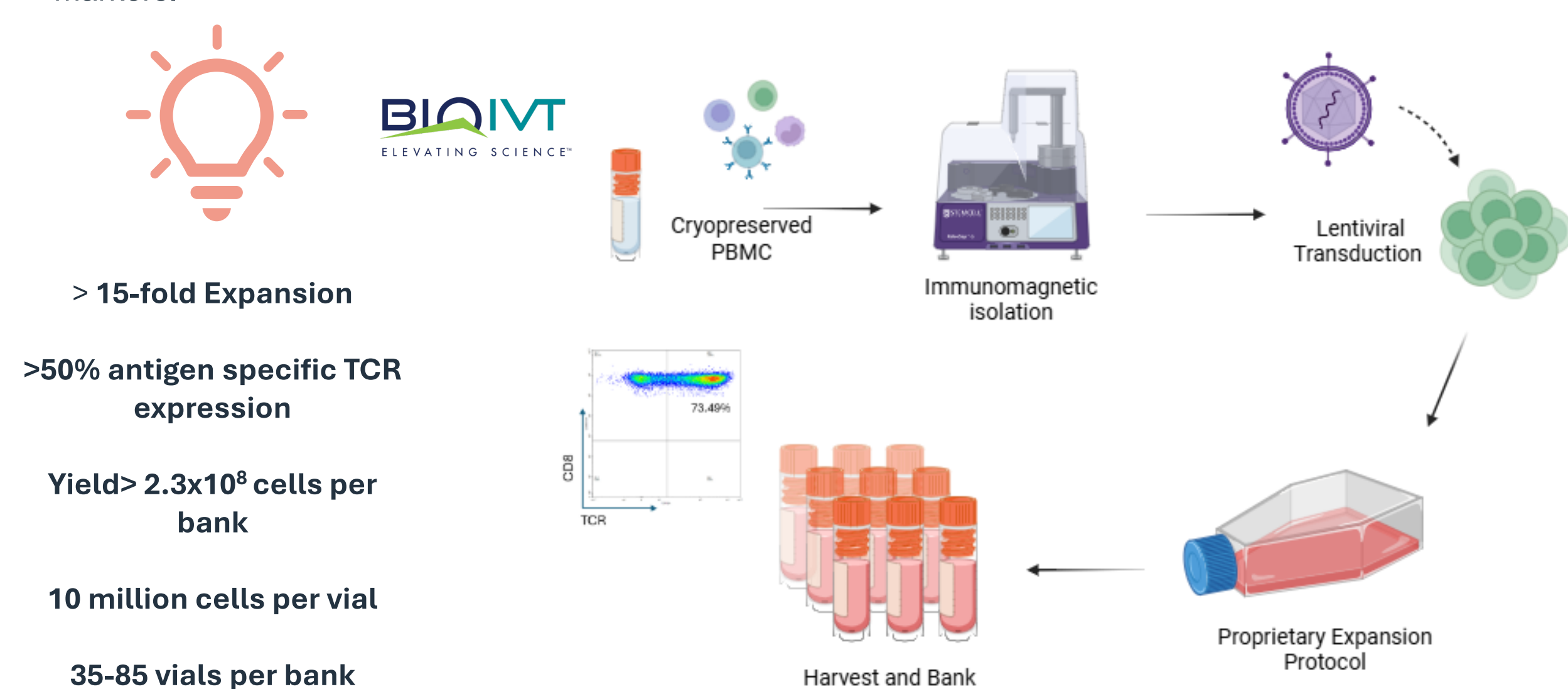


TCR-T Cells for Assessment of Immune Modulators

TCR-T cells were employed to evaluate the effects of immune modulators. T2 cells, which express HLA-A2 and can present specific antigens to cytolytic CD8+ T cells, were employed as the target cells. The T2 cells were pulsed with the antigen of interest and co-cultured with TCR-T cells in the presence of a concentration series of T cell inhibitors, enabling functional assessment of immune modulation.

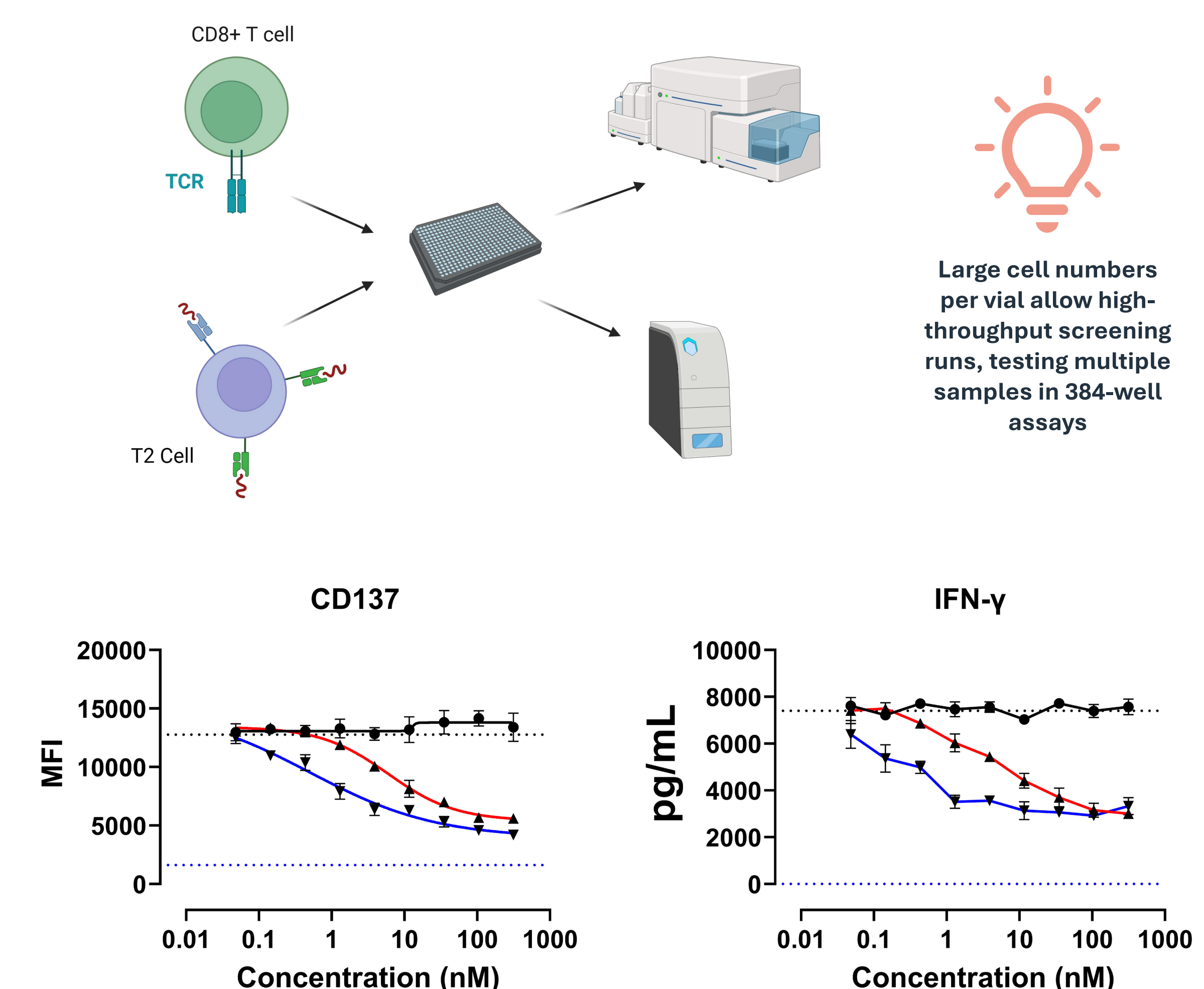


Functional Assessment of TCR-T Cells

MeWo cells are fibroblast-like cells isolated from a patient with malignant melanoma and have been reported to express the melanoma-associated antigen, MART-1 (Melan A). These cells were employed as target cells and co-cultured with CD8+ T cells engineered to express a transgenic MART-1-specific TCR at two E:T ratios (1:1 and 5:1), enabling antigen specific recognition and functional assessment of the TCR-T cells.

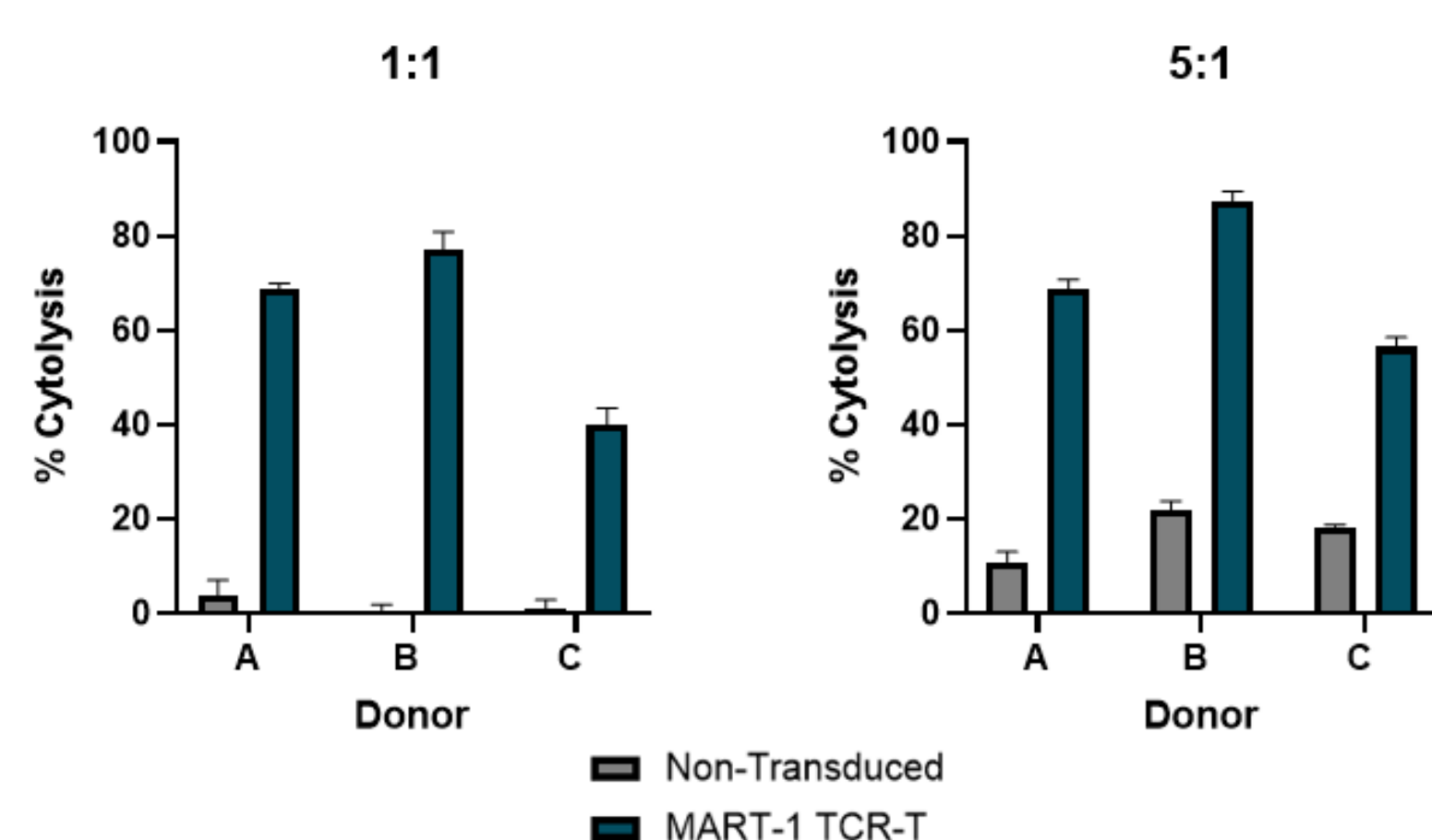


T2 cells were pulsed with either a MART-1 peptide, or an alternative non-specific peptide to evaluate antigen cross-reactivity. These peptide-loaded T2 cells were co-cultured with CD8+ T cells expressing a transgenic TCR specific for MART-1. T cell activation was assessed by measuring CD137 expression on CD8+ TCR+ cells, as well as assessing IFN-γ production.

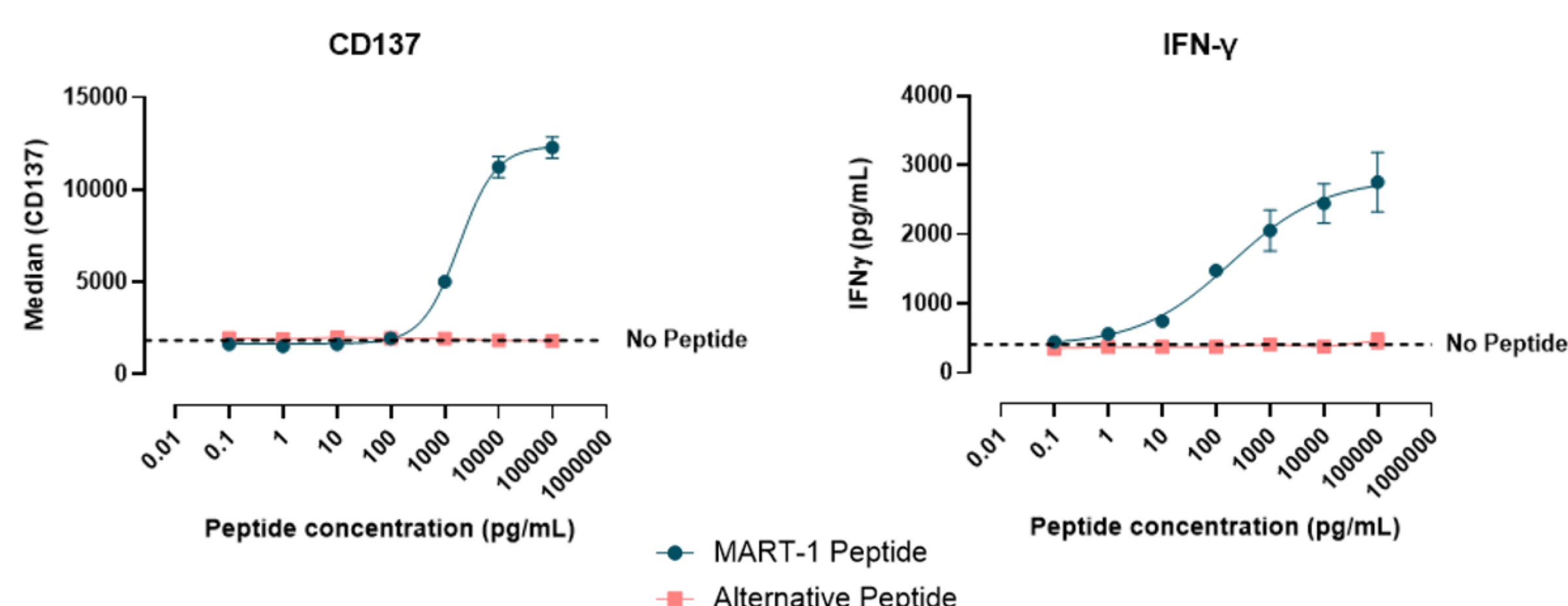

Endogenous Antigen-Specific T Cells

An alternative strategy is to generate antigen-specific T cells by expanding CD8+ T cells that endogenously express a TCR specific to the target antigen. These cells offer enhanced biological relevance by maintaining native TCR expression, without the need to introduce a second TCR.

Naïve CD8+ T cells were isolated from cryopreserved PBMC and expanded following a proprietary protocol. Functional activity was assessed using HLA-A1+ K562 cells pulsed with CMV peptide as antigen-presenting targets. Co-culture with CMV-specific endogenous antigen-specific T cells enabled evaluation of antigen-specific response.

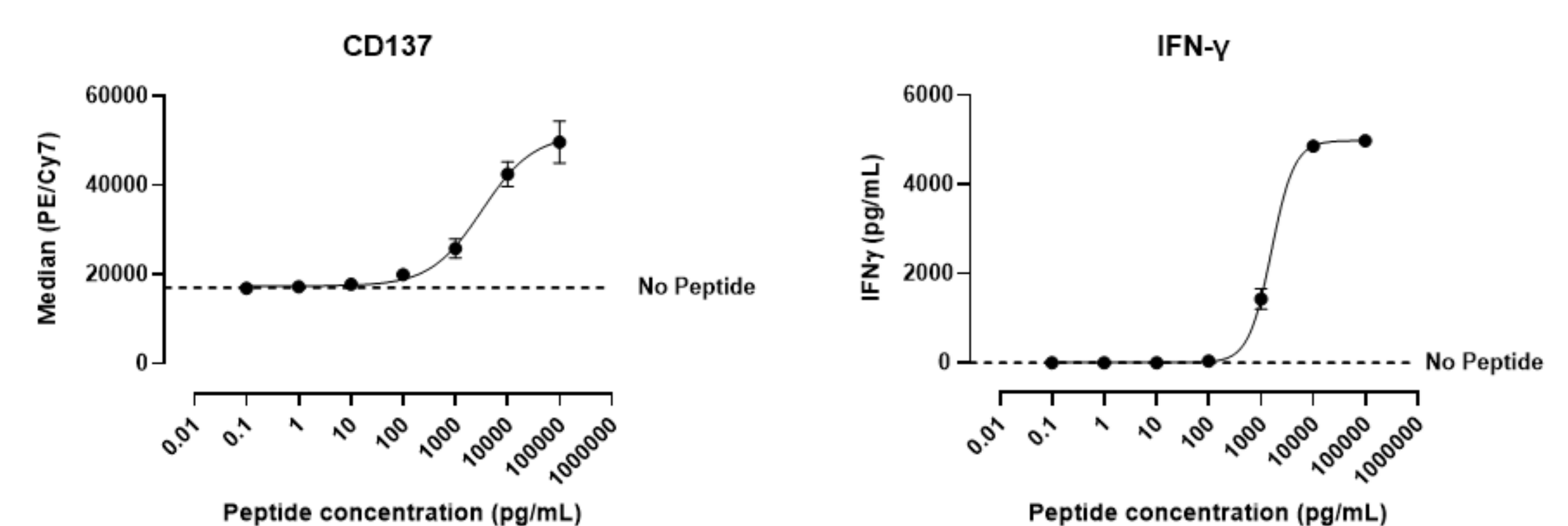

Summary

We have created 'Thaw & Go' banks of TCR-T cells (from PBMC supplied by BioIVT, LLC), ideal for high-throughput evaluation of immune-modulating therapies. These well-characterised cells exhibit specificity to their target antigen and can be assessed through a range of functional assays. Moreover, we have established an effective approach to expand endogenous antigen-specific T cells, providing a biologically relevant alternative. Finally, we are currently producing further cell banks, including Tregs and $\gamma\delta$ T cells, to offer valuable resources for a wider range of applications and to support faster candidate selection and advance drug discovery efforts.

Scan the code

Download the digital version of this research poster and take all the latest advancements and data home in your pocket.

